



253D Operation Manual



1. Dimension

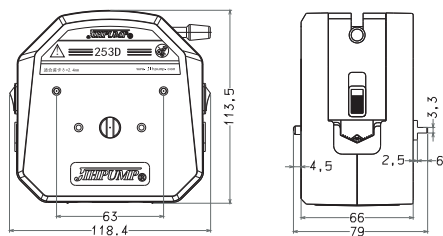

2. Flow rate

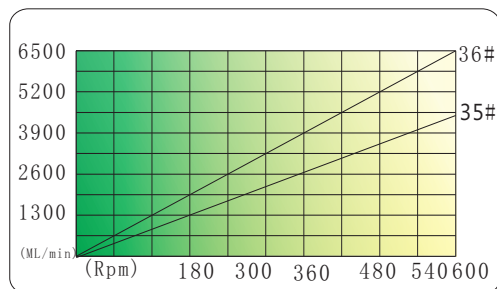

Model#	Tube	Housing	Roller	Pump case	Max flow	Weight
WT-600CA/253D	35# 36#	PPS& 30 GF	304	Plastics	6500mL/min	5.0Kg
WT-600CAS/253D				Stainless steel		6.35Kg
WT-600CL/253D				Cast Al alloy		6.4Kg
253D-ZL						2.8Kg



WT-600CA/253D



WT-600CAS/253D




WT-600EL/253D




253D-ZL


3. Install & use

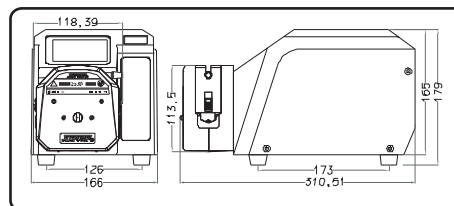


WT-600CA/253D : large LCD screen, basic and advanced data showed simultaneously, brushless motor drive, plastic housing.

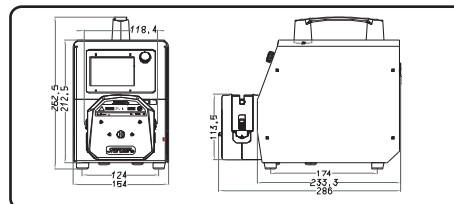

WT-600CAS/253D : large LCD screen, basic and advanced data showed simultaneously, brushless motor drive, stainless housing.

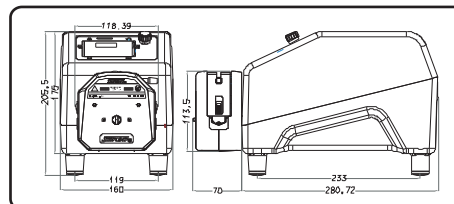

WT-600EL/253D : LED screen, brushless motor drive, cast Al alloy overall case, adapt to various bad working environment.

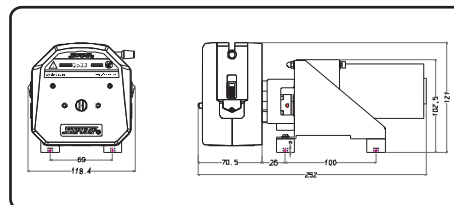

253D-ZL : DC motor drive, easy load type, use individually or combined use with instrument or equipment, with speed adjusted or constant.

4. Assemble tube



Ccw rotate lift pole



Get out press board



Assemble tube, press down board



Cw rotate lift pole, press tightly